

Gyrolab® Assays

Kineret® (anakinra) PK Assay


INTRODUCTION

Kineret (anakinra) is an anti-inflammatory biopharmaceutical drug used to treat rheumatoid arthritis and similar inflammatory diseases. It is a recombinant modified version of the human interleukin 1 receptor antagonist (IL-1 RA) protein and works by blocking the action of IL-1.

We have developed a three-step bridging Gyrolab PK Assay to determine Kineret in human serum samples. An MRD of 1:10 gives a broad analytical range with an approximate LLOQ of 50 ng/mL, and ULOQ of 100 000 ng/mL in neat serum. Use of this protocol on Gyrolab systems will reduce time to market and increase productivity while maintaining quality requirements.

ASSAY DESIGN

The assay was set up as a three-step bridging assay with biotinylated polyclonal anti-IL-1 RA as a capture molecule and monoclonal anti-IL-1 RA labeled with Alexa Fluor® 647 as a detection molecule.

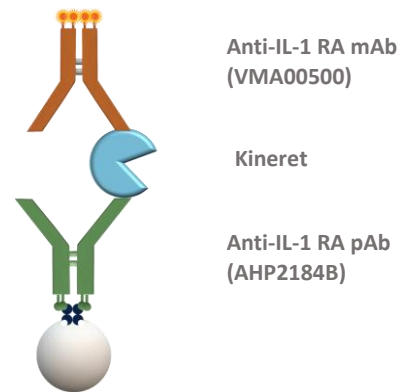

ASSAY PERFORMANCE

Dynamic range, accuracy and precision

A robust 3-log standard curve (Figure 1) was generated over three runs, achieving an assay range from 50 ng/mL to 100 000 ng/mL (Table 1). The inter-run precision (CV, Coefficient of Variation), established with QC samples over the assay range run in triplicate in three runs, was <20% (Table 2).

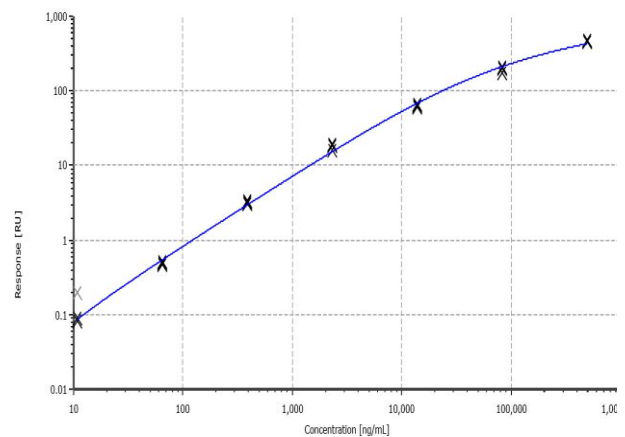


Figure 1 Standard curve in Rexasip H with 10% serum. Concentrations in neat serum

Table 1 Estimated Assay Range in neat serum, based on three runs

LLOQ (ng/mL)	ULOQ (ng/mL)
~ 50	~ 100 000

Table 2 Accuracy and precision data of QC samples in neat serum (n = number of runs)

QC	Expected Conc (ng/mL)	Average Measured Conc (ng/mL)	Inter Run CV (%; n=3)	Average Intra Run CV (%; n=3)	Average Total Error (%; n=3)
1	50	48	14	3.5	13
2	400	474	14	5.0	23
3	2 500	2 936	11	3.7	21
4	15 000	15 500	18	17	26
5	100 000	109 191	10	6.2	18

Selectivity

Selectivity was established by spiking 60 ng/mL of the drug in human serum samples. All samples measured <LLOQ when analyzed unspiked.

Table 3 Selectivity spiked samples

Sample	Expected Conc (ng/mL)	Average Measured Conc (ng/mL)	CV (%)	Average Bias (%)
1	60	58	7.6	-3.8
2	60	71	11	18
3	60	64	5.7	7.3
4	60	61	9.6	1.0
5	60	59	9.2	-1.6
6	60	62	5.9	2.5
7	60	60	4.2	0
8	60	64	2.6	7.0
9	60	58	5.3	-3.6
10	60	56	2.6	-6.5

MATERIALS AND METHODS

The assay was developed on a Gyrolab xP system using Gyrolab Bioaffy 1000 CD. The assay was set up using a 3-step Gyrolab method with two wash solutions (1000-3W-006-A) and a 1% PMT setting. The assay buffer was REXXIP H with 10% human serum. Anti-human IL-1 RA from BioRad (AHP2184B) was supplied biotinylated and used at a concentration of 700 nM, diluted with PBS-T.

The detection antibody labeled with Alexa Fluor 647 according to the Gyrolab standard protocol (Gyrolab User Guide), was monoclonal anti-IL-1 RA (BioRad, VMA00500), diluted to 35 nM in REXXIP F. The assay standard was the recombinant IL-1 RA anakinra from Swedish Orphan Biovitrum (Sobi). The standard was prepared in 10% human serum diluted in REXXIP H.

Summary table

Capture	700 nM biotinylated polyclonal anti-IL1 RA (AHP2184B, BioRad), in PBS-T
Detection	Alexa Fluor 647 labeled monoclonal anti-IL1 RA (VMA00500, BioRad) 35 nM in REXXIP F
Analyte	Kineret (Sobi) in REXXIP H with 10% human serum
CD-type	Bioaffy 1000 CD
Method	1000-3W-006-A
Wash buffer for needles	Wash buffer 1: PBS-T Wash buffer 2: Gyrolab Wash Buffer pH 11
PMT-setting	1%
Expected dynamic range	Approximately 50-100 000 ng/mL in neat human serum

Recommendations

When developing this assay for a specific drug development purpose, it is important to screen matrices and assess backgrounds, in particular for the specific disease matrices. Parameters, such as LLOQ should be validated in-house. Data given in this document should only be considered as guidance.

For additional support contact your local Field Application Support

Gyrolab and REXXIP are registered trademarks and Gyros, Gyrolab xPlore, Gyroplex, Bioaffy and Gyros logo are trademarks of Gyros Protein Technologies Group. All other trademarks are the property of their respective owners. Products and technologies from Gyros Protein Technologies are covered by one or more patents and/or proprietary intellectual property rights. All infringements are prohibited and will be prosecuted. Please contact Gyros Protein Technologies AB for further details. Products are for research use only. Not for use in diagnostic procedures. © All infringements are prohibited and will be prosecuted. Please contact Gyros Protein Technologies AB for further details. Products are for research use only. Not for use in diagnostic procedures. © Gyros Protein Technologies AB 2020. D0035600/A